

## Porta Wheels PSA-114

### INTRODUCTION:

PSA-114 Porta Wheels fit either the Gilson Porta-Screen® (PS-3 and PS-4) or the Gilson Sample Splitter (SP-1). Wheels provide mobility in the laboratory.

### ASSEMBLY:

Porta-Screens and Sample Splitters are supplied with mounting holes for fast installation of the Porta Wheels.

#### *To mount Porta Wheels on either machine:*

1. Insert cap screw through angle mount of wheel assembly.
2. Insert cap screw through predrilled mounting hole in machine, keeping the axle of wheel above mounting hole.
3. Install lock washer and nut, and tighten.



PSA-114



SP-1 shown with PSA-114



PS-4 shown with PSA-114

Rev: 06/20/2011